PRODUCT INFORMATION FOR USERS LEAFLET

1

STEEL-BRIGHT[®]MC

Sterile Pharmaceutical Clean Room Formula

Cleaner/Polish

| | | |
|--|--------------------------------------|------------------|
| Product Labeling | English | Page 2 |
| Le produit étiquète | Français(French) | Page 8 |
| Identificare del prodotto | Italiano(Italian) | Pagina 15 |
| Produkt-Beschriften | Deutsch(German) | Seite 22 |
| Etiquetado del producto | Español (Spanish) | Página 29 |
| 产品标签 | 中文 (普通话) Chinese (Mandarin) | 页36 |
| Märka för produkt | Svenska (Swedish) | Sida 43 |
| De Etikettering van het product | Nederlands (Dutch) | Pagina 50 |
| ملع عي جوتنم | [ك ي برأ] (Arabic) | صفحة 57 |
| Produkt maerke | Dansker (Danish) | Side 64 |

Veltek Associates, Inc.

15 Lee Boulevard • Malvern, PA 19355-1234

Tel: (610)644-8335 Fax: (610)644-8336 E-mail: vai@sterile.com PIUL STEEL-1002 EU/AP 7 JAN 2010

PRODUCT INFORMATION FOR USERS LEAFLET

2

STEEL-BRIGHT^{®MC}

Sterile Pharmaceutical Clean Room Formula

Read the label and leaflet before using

PRODUCT DESCRIPTION

STEEL-BRIGHT is an emulsion based professional strength stainless steel cleaner/polish. With STEEL-BRIGHT, stainless steel can be restored to a residue free finish. Brightens and polishes stainless steel. Removes spots created by water and oils. Will not rainbow or accumulate a heavy build up. Surfaces will remain cleaner longer because there is no residue film to attract soil. The gloss is renewed and will be retained. Product will not leave a powdery residue. This product can also be used on chrome, brass, copper and aluminum.

STEEL-BRIGHT Sterile Stainless Steel Cleaner addresses the use of a sterile cleaner/polish within the aseptic manufacturing area. The product is aseptically filtered, double bag packaged, and then gamma irradiated. The product is delivered each time with a lot specific Sterile Documentation Package. This innovation represents another significant breakthrough for clean room operation.

APPLICATIONS: For use on stainless steel or other smooth metals such as hand rails, light switch plates or shades, desks, stainless steel trim, wall or floor panels and doors, restaurant and cafeteria equipment. Can also be used on cars, trucks, motorcycles, engines, bicycles, boats, trailers, sporting equipment, guns, knives, jewelry, and many other applications.

- Removes spots and stains
- Brightens and Polishes
- Contains no acids or abrasives
- May be used on stainless steel, chrome, brass, copper & aluminum

General Product Information:

| | |
|--------------------------------|---|
| FILLING: | Product is filled in ISO 5 clean room Class 100 conditions. |
| FILTRATION: | Filtration designated per product is accomplished and integrity of the filter assured by prior and post bubble point testing |
| PRODUCT LOT NUMBER: | Product Fill Lot Number is placed on the bottom of each can, and on the exterior of the case. |
| IRRADIATION LOT NUMBER: | Irradiation Sterilization Lot Number appears on the exterior of the case. Dosage is Min: 25.0 kGy Max: 50.0 kGy |
| CAN: | Seamless Aluminum, cleaned to class 100, exterior wiped with IPA before packaging. |
| PACKAGING: | Product is double bagged per can in 4 mil polyethylene bags and placed into lined cases. The entire contents are irradiated. |
| STERILITY TESTING: | Once irradiation treatment has been accomplished, the product undergoes 14 day USP Sterility Test - Direct Transfer Method (20 samples/lot) |

Veltek Associates, Inc.

15 Lee Boulevard • Malvern, PA 19355-1234

Tel: (610)644-8335 Fax: (610)644-8336 E-mail: vai@sterile.com PIUL STEEL-1002 EU/AP 7 JAN 2010

PRODUCT INFORMATION FOR USERS LEAFLET

4

STEEL-BRIGHT[®]MC

PHYSICAL AND CHEMICAL HAZARDS

This product may be irritating to the eyes, skin, gastrointestinal tract and respiratory system. Contents under pressure. Do not puncture or incinerate container. Exposure to temperatures above 48.8°C (120° F) may cause bursting. Direct contact with aerosol propellant may cause frostbite or freezing of exposed skin. Avoid puncture, creasing or impacting side walls of aerosol container.

Skin Effects

This product may cause irritation to the skin.

Eye Effects

This product may cause irritation to the eyes.

Ingestión Effects

Ingestión of this product may result in central nervous system effects including headache, sleepiness, dizziness, slurred speech and blurred vision. Material, if swallowed, can be aspirated into lungs.

Inhalation Effects

May cause irritation to the nose and respiratory tract. Inhalation may cause central nervous system depression with symptoms such as weakness, dizziness, confusion and drowsiness.

PERSONAL PROTECTIVE EQUIPMENT:

Eyes/Face

Wear chemical goggles; face shield (if splashing is possible).

Skin

Chemical resistant gloves should be used at all times when handling this product.

Respiratory

If ventilation is not sufficient to effectively prevent buildup of vapors, appropriate NIOSH/MSHA respiratory protection must be provided

Personal Protective Equipment: General

Eye wash fountain and emergency showers are recommended.

Chemical Stability

Stable under recommended conditions of storage and handling.

Conditions to Avoid: Keep away from heat, ignition sources and incompatible materials.

Incompatibility: Acids, oxidizing materials, combustible materials, peroxides, chlorine, and reactive metals.

Hazardous Decomposition: Oxides of carbon.

Possibility of Hazardous Reactions: Will not occur

Veltek Associates, Inc.

15 Lee Boulevard • Malvern, PA 19355-1234

Tel: (610)644-8335 Fax: (610)644-8336 E-mail: vai@sterile.com PIUL STEEL-1002 EU/AP 7 JAN 2010

PRODUCT INFORMATION FOR USERS LEAFLET

5

STEEL-BRIGHT[®]MC

DIRECTIONS for USE

Shake can well before using. Hold can upright, spray a small amount to wet surface. Clean with one cloth and polish with another clean, dry cloth. DO NOT use on pots, pans or eating utensils. [Note: The use of this product on sterile surfaces in aseptic controlled areas should always be done by trained individuals using good aseptic technique and sterilized cloths or wipes. In non-sterile areas, clean non-sterile materials may be used for application and polishing.]

SAFETY HAZARD INFORMATION

For a complete summary of the chemical and physical safety and exposure hazards to users, see **Material Safety Data Sheet SB-01112**. A copy of the MSDS was shipped with this product. If an additional copy is needed contact VAI USA at **+610-644-8335 or FAX +610-644-8336** Veltek Customer Service or Website www.sterile.com (If an immediate copy is needed, outside business hours, call the Emergency Response Service listed below.)

(Emergency): For Spill/Exposure Emergency Response Service from Europe in English (and 23 European languages) telephone CARECHEM24 at +44 1235 239 670, for Arabic language +44 1235 239 671, for Chinese +86 10 5100 3039

EU Classification and Labelling according to Directive(s) 67/548/EEC and/or 1999/45/EC:

Symbol(s)

Xn Harmful Risk

Phrases

R65- Harmful: may cause lung damage if swallowed.

Safety Phrases

S23 - Do not breathe gas/fumes/vapour/spray.

S24 - Avoid contact with skin.

S62- If swallowed, do not induce vomiting. Seek medical advice immediately and show the container or label.

Substance/Preparation Classification according to CLP (EC) Regulation No. 1272/2008:

Aspiration Hazard; Category 1

Acute Toxicity (dermal); Category 5 (90% unknown)

Acute Toxicity (inhalation); Category 5 (90% unknown)

Acute Aquatic Toxicity; Category 3

Veltek Associates, Inc.

15 Lee Boulevard • Malvern, PA 19355-1234

Tel: (610)644-8335 Fax: (610)644-8336 E-mail: vai@sterile.com PIUL STEEL-1002 EU/AP 7 JAN 2010

PRODUCT INFORMATION FOR USERS LEAFLET

6

STEEL-BRIGHT^{®MC}

Label Elements

Signal Word

DANGER

Hazard Statement(s)

May be fatal if swallowed and enters airways.

May be harmful in contact with skin.

May be harmful if inhaled.

Harmful to aquatic life.

Precautionary Statement(s)

Prevention

Do not spray on an open flame or other ignition source. Pressurized container: Do not pierce or burn, even after use. Avoid release to the environment.

Response

Call a POISON CENTER or doctor/physician if you feel unwell.

IF INHALED: Call a POISON CENTER or doctor/physician if you feel unwell.

IF SWALLOWED: Immediately call a POISON CENTER or doctor/physician. Do NOT induce vomiting.

Repeated Dose Toxicity

Repeated or prolonged exposure may cause skin disorders including eczematous or contact dermatitis, folliculitis, oil granuloma, and melanosis.

STORAGE AND DISPOSAL

Do not contaminate water, food, or feed by storage and disposal. Keep out of reach of children.

Storage

Store in the original container in a dry cool place. Protect from sunlight. Do not expose to temperatures exceeding 50°C (122°F).

General Product Information

In the US A, Wastes must be tested using methods described in 40 CFR Part 261 to determine if it meets applicable definitions of hazardous wastes.

Component Waste Numbers

No USA EPA Waste Numbers are applicable for this product's components.

Disposal Instructions

Dispose of contents/container in accordance with application regulations.

Veltek Associates, Inc.

15 Lee Boulevard • Malvern, PA 19355-1234

Tel: (610)644-8335 Fax: (610)644-8336 E-mail: vai@sterile.com PIUL STEEL-1002 EU/AP 7 JAN 2010

PRODUCT INFORMATION FOR USERS LEAFLET

7

STEEL-BRIGHT[®]MC

Container Disposal

Do not reuse or try to refill this container. Do not puncture container. Do not incinerate container. Offer for recycling, or dispose of in a sanitary landfill.

If you have any questions, your local VAI Technical Support Representative can assist you in assuring all considerations are addressed prior to use. Please contact VAI USA at **+610-644-8335, FAX +610-644-8336** Veltek Customer Service for the name and number of your local VAI Sales or Technical Services representative or www.sterile.com.

LOT SPECIFIC STERILE DOCUMENTATION (delivered with each shipment)

- Certificate of Analysis
- Certificate of Sterility
- Certificate of Irradiation

(Please contact VAI for a sample of this documentation)

Available Technical Data Supplements (Upon Request)

LUBE-1001: Sterile Lubricant and Cleaner Guide

Please contact VAI USA at **+610-644-8335, FAX +610-644-8336** Veltek Customer Service for the name and number of your local VAI Sales or Technical Services representative for additional copies.

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